

Synchronize your controller to your Standard

This procedure will synchronize your controller with your Standard.

1. Turn the controller ON and let it pass the SELF TEST.
2. Press and hold the ENTER (right most key) and the UP-ARROW key TOGETHER for 1 second. (Out 1 should show on the bottom of the display).
3. Scroll down using the INDEX key (left most key) until you reach InPC.
4. Change the value positive or negative to match the DIFFERENCE in the temperature using the UP or DOWN ARROW keys. (It can be adjusted + or - up to 500 degrees.)
5. Press the ENTER key (right most key). (Display should blink.)
6. Turn the unit OFF and back ON.

This completes the Synchronization process.

If the Controller number does not match your Standard, you can go back and repeat Steps 1 thru 6 until it does.